



BIOTECH AB

WWW.BIOTECH-SWEDEN.COM

Application Data


No. 1013A

ChromaNik
ChromaNik Technologies Inc.

バルサルタンと関連物質の分離

Separation of Valsartan and Related Substances


Sunniest C18 3 μ m 150 x 4.6 mm i.d.

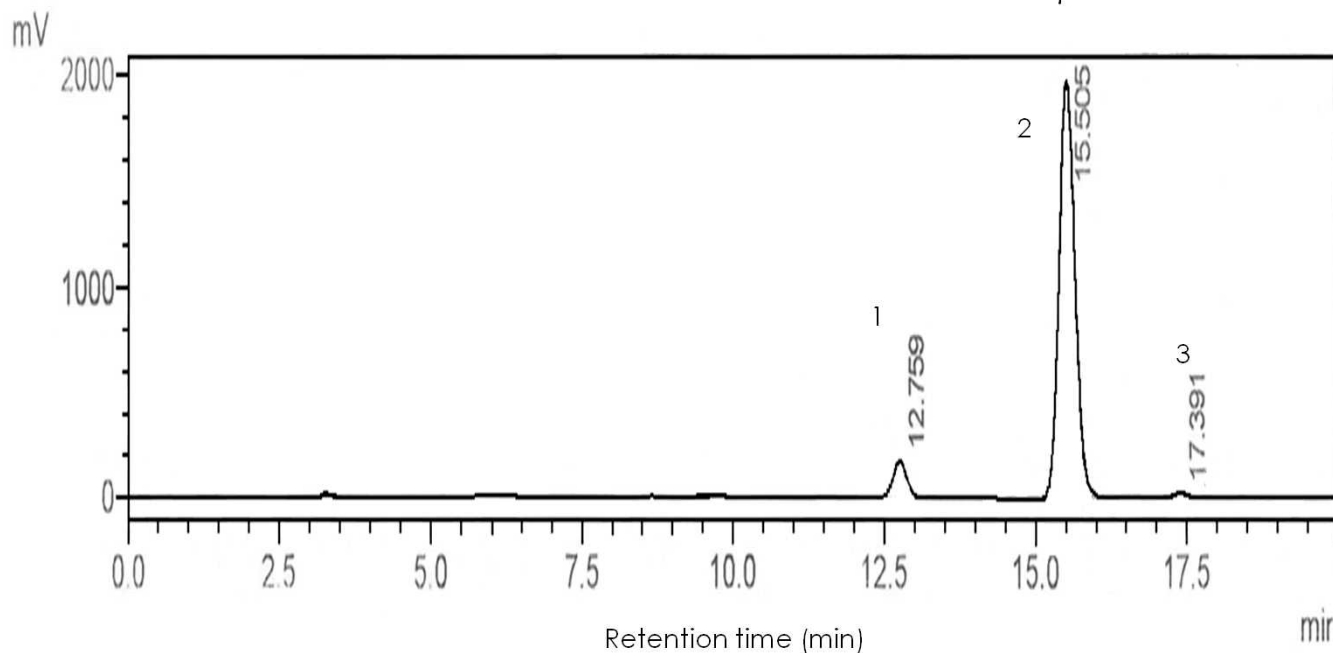

1 Unknown

2

3 Unknown

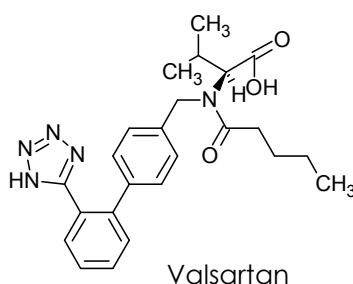

Column: Sunniest C18, 3 μ m 150 x 4.6 mm

Mobile phase: Acetonitrile/Water/Glacial Acetic Acid = 500/500/1

Flow rate: 0.4 mL/min

Temperature: 30 °C

Detection: UV@225nm

Test probe : Valsartan 0.5mg/mL in mobile phase

Injection volume: 10 μ L

Biotech AB (Sweden) Email: info@biotech.se

Biotech USA Email: info@biotechusa.us

Biotech India Pvt Ltd Email: tech@biotechltd.in

ChromaNik Technologies Inc. (Japan) Email: info@chromanik.co.jp

No. 1013A