


## Pharmaceuticals


## Depressants

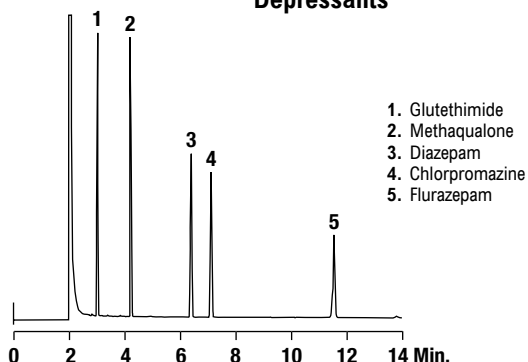

**Column:** Heliflex® AT™-5ms, 30m x 0.25mm x 0.25µm (Part No. 15807)  
**Temp:** 270°C  
**Carrier Gas:** Helium at 0.65mL/min (24cm/sec)  
**Detector:** FID

CHROM 2560


## Anticonvulsants

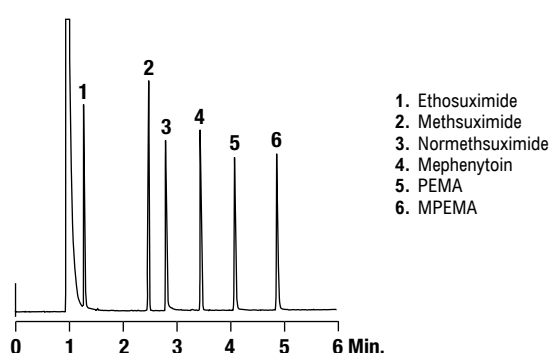

**Column:** Heliflex® AT™-5ms, 15m x 0.25mm x 0.25µm (Part No. 15801)  
**Temp:** 190°C (1 min hold) to 235°C at 10°C/min  
**Carrier Gas:** Helium at 0.7mL/min (26cm/sec)  
**Detector:** FID

CHROM 2568


## Antidepressants

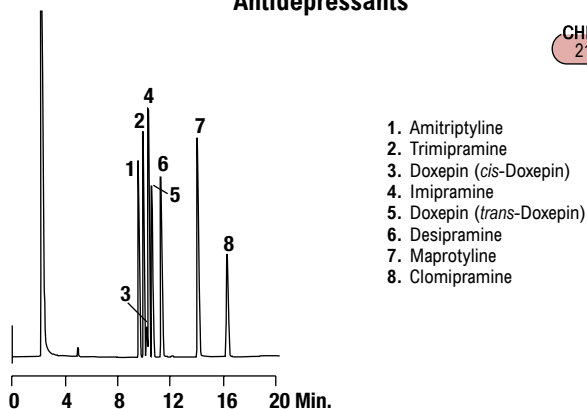

**Column:** Econo-Cap™ AT™-20, 30m x 0.32mm x 0.25µm (Part No. 19677)  
**Temp:** 235°C  
**Carrier Gas:** Helium at 22 cm/sec  
**Detector:** FID

CHROM 2187


## Antihistamines

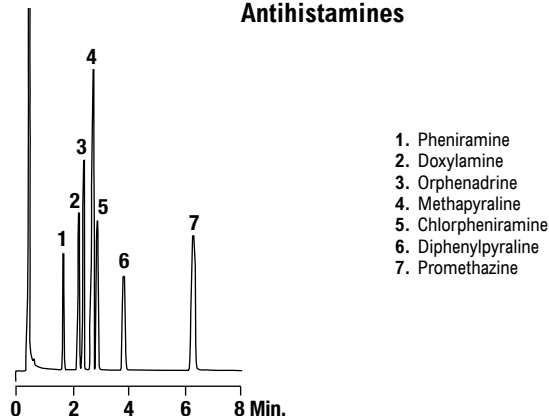

**Column:** Econo-Cap™ EC™-1, 15m x 0.53mm x 1.20µm (Part No. 19650)  
**Temp:** 210°C  
**Carrier Gas:** Helium at 8.5mL/min  
**Detector:** FID

CHROM 1005


## Antihistamines

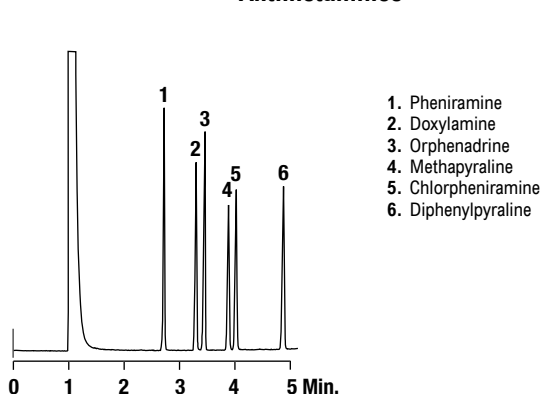

**Column:** Heliflex® AT™-5ms, 15m x 0.25mm x 0.25µm (Part No. 15801)  
**Temp:** 210°C (1 min hold) to 230°C at 5°C/min  
**Carrier Gas:** Helium at 0.7mL/min (25cm/sec)  
**Detector:** FID

CHROM 2567


## Hypnotics

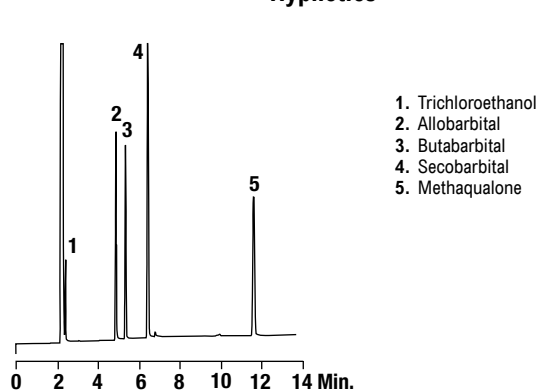

**Column:** Econo-Cap™ EC™-1, 30m x 0.32mm x 0.25µm (Part No. 19651)  
**Temp:** 160°C to 220°C at 10°C/min  
**Carrier Gas:** Helium at 1.08mL/min  
**Detector:** FID


CHROM 1009

## Fast GC Drug Analysis

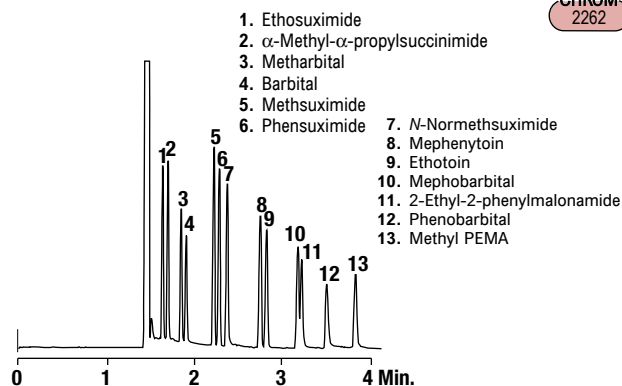CHROM  
2262


Column: Heliflex® AT™-5, 20m x 0.10mm x 0.10 $\mu$ m  
(Part No. 14454)  
Temp: 240°C  
Carrier Gas: Helium at 0.1mL/min  
Detector: FID

## Fluoxetine and Metabolite

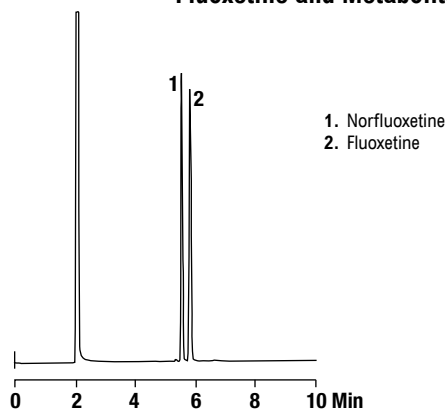CHROM  
1786


Column: Heliflex® AT™-1, 30m x 0.32mm x 0.25 $\mu$ m  
(Part No. 13640)  
Temp: 200°C  
Carrier Gas: Helium at 1.06mL/min  
Detector: FID

## Sedatives

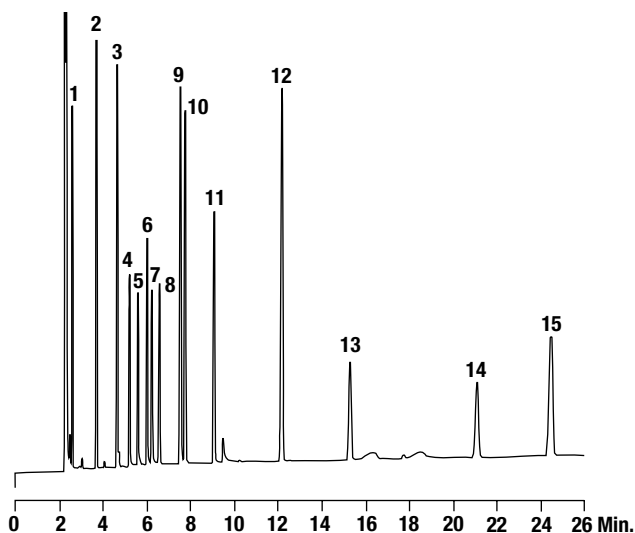CHROM  
1003

Column: Econo-Cap™ EC™-1, 30m x 0.32mm x 0.25 $\mu$ m  
(Part No. 19651)  
Temp: 160°C to 220°C at 10°C/min  
Carrier Gas: Helium at 1.08mL/min  
Detector: FID

## Nitrazepam and ANB Metabolite

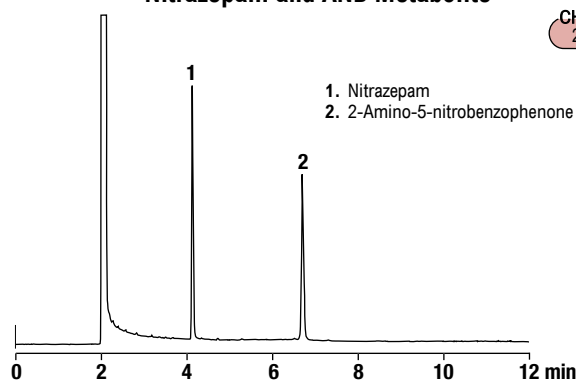CHROM  
2592

Column: Heliflex® AT™-1ms, 30m x 0.25mm x 0.25 $\mu$ m  
(Part No. 15897)  
Temp: 280°C  
Carrier Gas: Helium at 0.7mL/min (25cm/sec)  
Detector: FID

## more applications

View our complete searchable chromatogram database at [www.discoverysciences.com/chromdb/](http://www.discoverysciences.com/chromdb/)