# Purification of Conjugated Quercetin with Rutinose Using Reveleris<sup>®</sup> Flash System is Fast and Productive

Reveleris® iES Flash Chromatography System

#### Introduction

Flavanol glycoside Rutin has been used in dietary supplements as Vitamin P for therapeutic treatments.<sup>1</sup> Besides its antioxidant property, this bioflavonoid works with vitamin C and helps in reducing pain and intraocular pressure. It has been synthesized from the conjugation of carbohydrates such as rhamnose and glucose with the flavonoid guercetin.


# Experimental

Run Conditions
Cartridge: Reveleris® amino 12g
Flow rate: 36 mL/min
Equilibration: 10.0 min
Run length: 9.0 min
UV1 wavelength: 265 nm
UV2 wavelength: 220 nm
ELSD carrier: Iso-proponal
Solvent A: Water
Solvent B: Acetonitrile with 0.5% TFA

<b>Gradient Method</b>		
Step	Time (min.)	%B
1	0	99
2	1	99
3	3	70
4	5	70



## **Results and Discussion**

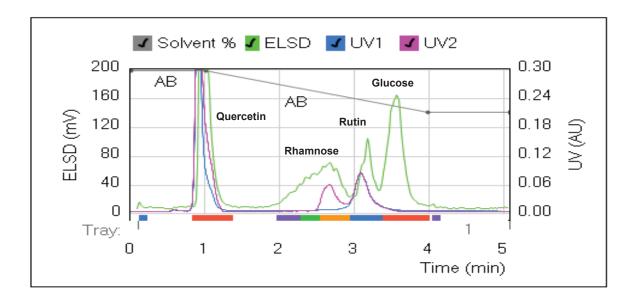

Rutin

### Conclusion

Using the Reveleris® amino cartridge and the ELSD detector of the RevealX™ technology, sugar impurities from rutin can be isolated and purified. With other lower capacity cartridges, closely eluting peaks may co-elute

resulting in poor recovery during isolation. As amino phase media is ideal for separating sugar compounds, all the components in the reaction mixture are separated using trifluoroacetic acid as the solvent modifier with speed.

#### References

1. Medicinal natural products; a biosynthetic approach, 3rd edition; Dewick, P.; John Wiley & Sons, Inc., Hoboken, New Jersey, 2009.

Americas Deerfield, Illinois Tel: +1 847.948.8600 discoverysciences@ grace.com

Europe Lokeren, Belgium Tel: +32 (0)9.340.65.65 discoverysciences.EU@ grace.com India Pune Tel: +91 20.6644.9900 pune@grace.com China Shanghai Tel: +86 21.5467.4678 dsbiz.asia@grace.com Australia Rowville Tel: +61 3.9237.6100 discoverysciences.AU@ grace.com

GRACE®, GRACE DAVISON® are trademarks, registered in the United States and/or other countries, of W. R. Grace & Co.-Conn. REVELERIS® is a trademark, registered in the United States and/or other countries, of Alltech Associates, Inc.

REVEALX™ is a trademark of Alltech Associates, Inc. This trademark ist has been compiled using available published information as of the publication date of this brochure and may not accurately reflect current trademark ownership or status.

Grace Daviscon Discovery Sciences is a product group of W. R. Grace & Co.-Conn., which now includes all product lines formerly sold under the Alltech Associates, Inc. is a wholly owned subsidiary of W. R. Grace & Co.-Conn.

The information presented herein is derived from our testing and experience. It is offered for your consideration and verification. Since operating conditions vary significantly, and are not under our control, we disclaim all warranties on the results that may be obtained from the use of our products. W. R. Grace & Co.-Conn. and its subsidiaries on to be held responsible for any damage or injury occurring as a result of improper installation or use of its products. Grace reserves the right to change prices and/or specifications without prior notification. © Copyright 2011 W. R. Grace & Co.-Conn. All rights reserved.

GRACE