

RevealX™ Technology Can be Used in the Impurity Isolation of Valproic Acid from Cyclodextrin During Encapsulation

Reveleris® iES Flash Chromatography System

Introduction

Valproic acid is one of the major antiepileptic drugs used in the treatment of different kinds of epilepsy.¹ The drug has shown to be effective in treating cancer due to its antitumor properties. Although water is usually used as the reaction medium for enzymes, organic solvents or

cyclodextrins have been added to improve the solubility of poorly soluble compounds. To increase its water solubility and bioavailability, the drug can be linked chemically by physical entrapment with polymers such as cyclodextrin or to polymeric carriers such as dextran.

Experimental

Run Conditions

Cartridge: Reveleris® C18 12g

Flow rate: 36 mL/min

Equilibration: 3.0 min

UV1 wavelength: 220 nm

UV2 wavelength: 254 nm

ELSD carrier: Iso-proponal

Solvent A: Water

Solvent B: Acetonitrile

Gradient Method

Step	Time (min.)	%B
1	0	30
2	7	80

Run Conditions

Cartridge: Reveleris® amino 12g

Flow rate: 36 mL/min

Equilibration: 6.0 min

Run length: 12.0 min

UV1 wavelength: 212 nm

UV2 wavelength: 254 nm

ELSD carrier: Iso-proponal

Solvent A: Water with 0.1% TFA

Solvent B: Acetonitrile

Gradient Method

Step	Time (min.)	%B
1	0	90
2	1	90
3	4	70
4	7	70



Results and Discussion



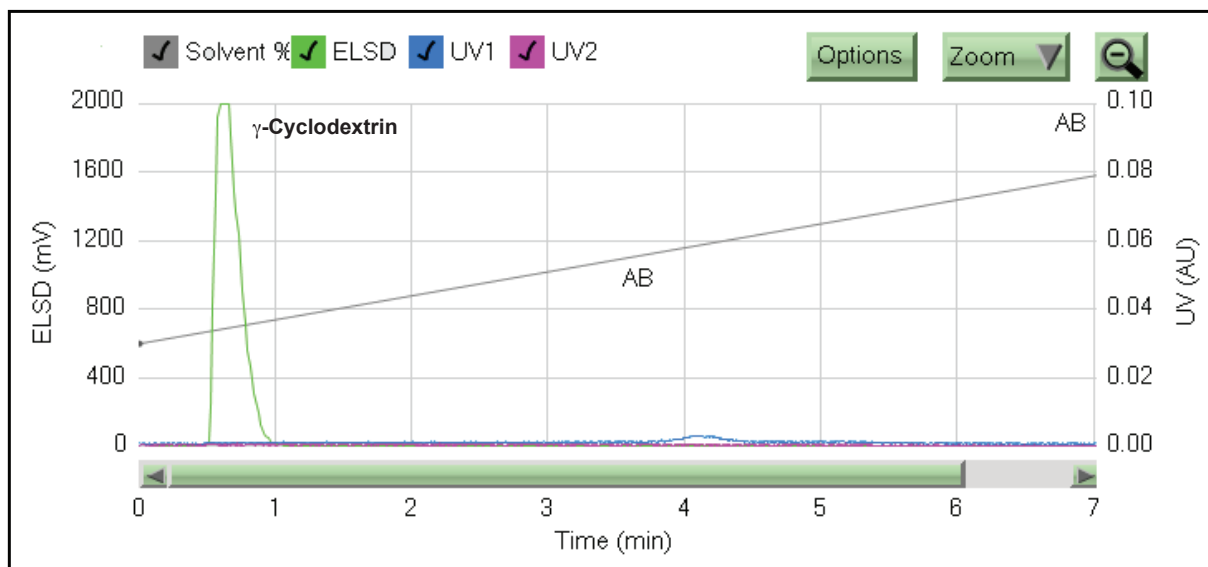


Figure 1: In this application, the separation of valproic acid from its entrapment-carrier, γ -cyclodextrin, has been shown using the Reveleris® flash cartridge with the RevealX™ detection technology. Using a linear gradient with acetonitrile and water mobile phase, cyclodextrin is poorly retained on a reversed phase cartridge.

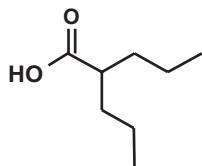

Valproic acid

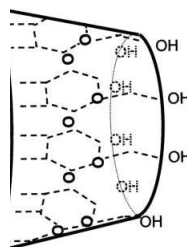

γ -Cyclodextrin



Results and Discussion

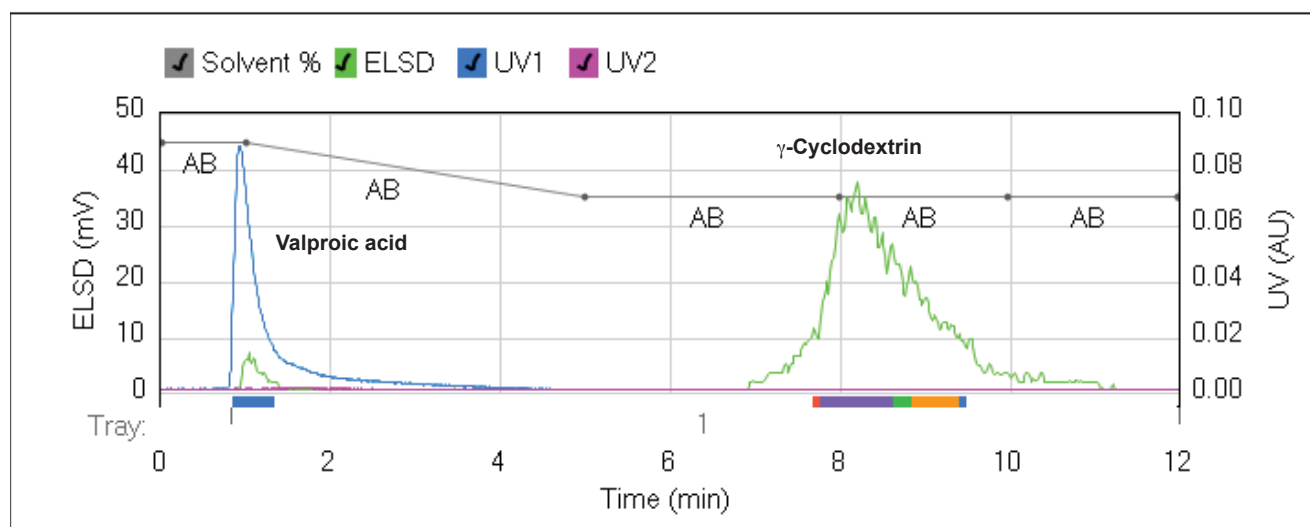


Figure 2: Compared to a reversed phase cartridge separation as shown where the retention is based on the hydrophobic interaction, the water-soluble cyclodextrin has been retained and separated from the free valproic acid using an amino phase media cartridge.

Conclusion

Polymer nanoparticles allow proper drug delivery and targeting with a wide range of properties under controlled conditions.² As polymer technology plays a key role in overcoming drug delivery challenges for those undeliverable molecules, nanospheres and nanocapsules designed as drug carriers can be synthesized with drugs linked through encapsulation or conjugation for better

target delivery. There needs to be a better purification technique for isolation of these compounds when the drug is poorly soluble in water while its carrier is a polar molecule. Hence such a method helps medicinal chemists isolate water based encapsulated drug from its impurity either as free or degraded and unbound compounds from the reaction mixture.

References

1. Praveen, B., Shrivastava, P., Shrivastava, S. K.; In-Vitro release and pharmacological study of synthesized valproic acid-dextran conjugate; *Acta Pharmaceutica Scientia*, (2009), 51, pp. 169 – 176.
2. Vauthier, C. and Bouchemal, K.; Methods for the preparation and manufacture of polymeric nanoparticles; *Pharmaceutical Research*, (2008), 26, No. 5, pp. 1025 – 1058.

Americas
Deerfield, Illinois
Tel: +1 847.948.8600
discoverysciences@
grace.com

Europe
Lokeren, Belgium
Tel: +32 (0)9.340.65.65
discoverysciences.EU@
grace.com

India
Pune
Tel: +91 20.6644.9900
pune@grace.com

China
Shanghai
Tel: +86 21.5467.4678
dsbiz.asia@grace.com

Australia
Rowville
Tel: +61 3.9237.6100
discoverysciences.AU@
grace.com

GRACE®, GRACE DAVISON® are trademarks, registered in the United States and/or other countries, of W. R. Grace & Co.-Conn. REVELERIS® is a trademark, registered in the United States and/or other countries, of Alltech Associates, Inc. REVEALX™ is a trademark of Alltech Associates, Inc. This trademark list has been compiled using available published information as of the publication date of this brochure and may not accurately reflect current trademark ownership or status. Grace Davison Discovery Sciences is a product group of W. R. Grace & Co.-Conn., which now includes all product lines formerly sold under the Alltech brand. Alltech Associates, Inc. is a wholly owned subsidiary of W. R. Grace & Co.-Conn. The information presented herein is derived from our testing and experience. It is offered for your consideration and verification. Since operating conditions vary significantly, and are not under our control, we disclaim all warranties on the results that may be obtained from the use of our products. W. R. Grace & Co.-Conn. and its subsidiaries can not be held responsible for any damage or injury occurring as a result of improper installation or use of its products. Grace reserves the right to change prices and/or specifications without prior notification. © Copyright 2011 W. R. Grace & Co.-Conn. All rights reserved.